



# LED Transilluminator, Triple wavelength

---

**Cat. No. BT-509**

This LED transilluminator is suitable for research use only.

It must be used by specialized personnel that know the health risks associated with blue light irradiation and with the reagents that are normally used with this instrument. Human naked eyes with long time irradiation of red, green, or blue light may have the probability of suffering from retina diseases. The amber filter provides blue light protection. The red filter provides green light protection. However, it does not guarantee complete protection. It is designed to shield the person working at the LED transilluminator only.

### **WARRANTY**

The LED transilluminator is warranted against defects in materials and workmanship for 1 year. If any defects occur in the instrument or accessories during this warranty period, BT Lab Systems will repair or replace the defective parts at its discretion without charge. The following defects, however, are specifically excluded:

1. Defects caused by improper operation.
2. Repair or modification done by anyone other than BT Lab Systems or an authorized agent.
3. Damage caused by substituting alternative parts.
4. Use of fittings or spare parts supplied by anyone other than BT Lab Systems.
5. Damage caused by accident or misuse.
6. Damage caused by disaster.
7. Corrosion caused by improper solvent or sample.

For any inquiry or request for repair service, contact BT Lab Systems. Inform BT Lab Systems of the model and serial number of your instrument.

### **REGULATORY NOTICE**

**IMPORTANT:** This BT Lab Systems instrument is designed and certified to meet safety standards and EMC regulations. Certified products are safe to use when operated in accordance with the instruction manual. This instrument should not be modified or altered in any way. Alteration of this instrument will:

1. Void the manufacturer's warranty
2. Void the safety and EMC certification
3. Create a potential safety hazard

BT Lab Systems is not responsible for any injury or damage caused by the use of this instrument for purposes other than those for which it is intended, or by modifications of the instrument not performed by BT Lab Systems or an authorized agent.

## IMPORTANT NOTICE

Please, read the installation instruction carefully before installing the LED transilluminator. This instrument is intended for clinical and research laboratory use with gel activation and it must be operated only by specialized personnel aware of the potential risks associated with the chemical and biological agents normally used with this unit. This instrument is meant for use only by specialized personnel that know the health risks associated with red, green, and blue light radiation and with reagents that are normally used with this instrument.

## SPECIFICATIONS

- Dimensions (mm): 280D x 340W x 80H
- Viewing surface (mm): 160D x 200W
- Wavelength (nm): 470, 530, 630
- LED arrangements: delta matrix for two-side illumination
- LED lifetime: 50,000 hours
- Power: 24V-1.25 A power adaptor
- Weight (Kg): 5

## INSTALLATION

Carefully unpack the transilluminator and the shield as follow:

1. First remove the bubble material at the top.
2. Remove the transilluminator from the two bubble material shells and place it on a stable, horizontal surface.
3. Remove the plastic protection film from the transparent glass viewing surface.

### ***Stand alone installation***

1. The instrument must be placed on a bench leaving at least 10 cm of space all around in order to avoid any obstacle that may reduce the ventilation.
2. Connect the instrument to the power using the annexed cable. The power font must be able to deliver at least 250 VA with a voltage between 100 and 240 Vac. The plug must have a ground connection. The adaptor supplies 24 V DC voltage.

### ***Installation with the Standard Documentation System:***

This instrument has been designed to work with the Standard Documentation System. In this case follow the instructions included in the Standard Documentation manual or any other instrument that will be released in the future.

## GENERAL PRECAUTIONS

- Plug the transilluminator on an electric line with ground connection.
- The transilluminator is equipped with thermal protection to prevent overheating.
- Do not pour liquids directly on the transilluminator.
- Do not block the aeration slits.
- Switch off the instrument immediately after its use.
- Position the transilluminator to prevent harm to nearby operators.
- When using the transilluminator with samples stained with specific dyes, decontaminate the transilluminator surface with bleach. Denatured alcohol can be also used. Always wear disposable gloves.

## USING THE TRANSILLUMINATOR

Place gel/sample on the transparent glass support area. It is recommended that gloves be worn to prevent skin contact with gel and staining agents. Press the ON/OFF switch to ON. The LEDs within the unit will begin glowing beneath the glass support. Dial the RGB knob to the appropriate light color that is intended to use. After viewing the sample, turn the transilluminator off via the ON/OFF switch and deplug the power adapter.

## GENERAL APPEARANCE OF THE LED TRANSILLUMINATOR



Adaptor input: 24V DC

Power AC input: 100~240 V; 50/60 Hz power source

## TECHNICAL SUPPORT

BT Lab Systems offers technical support for all of its products. If you have any questions about the product's use or, operation, please contact BT Lab Systems at the following info.

E-Mail: [info@BTLabSystems.com](mailto:info@BTLabSystems.com)